

Application Note

EVUSEP

Unlocking the plasma proteome with Evotip Pure based, standardized and scalable workflows

Highlights

- Fully automated, end-to-end workflow for neat and deep plasma profiling
- Standardized and cost-efficient strategy to monitor disease at scale

1. Introduction

Plasma stands out as one of the most readily accessible biofluids, with an estimated presence of over 10,000 distinct proteins. A significant number of these are released from organs and tissues. The ability to measure such proteins in a robust manner holds tremendous potential for advancing next-generation multiplexed biomarker discovery. However, methods must be able to deal with dynamic range of 10-12 orders of magnitude, while providing sufficient throughput in a robust and cost-effective manner. While a standardized approach is needed for wide utilization in the field, complementary strategies are also relevant to apply, depending on the use case. Here, we describe a fully automated workflow on the Opentrons OT-2 liquid handling robot (OT-2) for neat plasma preparation based on Protein Aggregation Capture (PAC)¹, and an approach for

deeper profiling of a sub-proteome based on the Mag-Net method². The workflows are completely hands-off, incorporating all steps from raw plasma to peptides loaded on Evotips. This is driven by efficient protocols based on hyper-porous magnetic beads directly followed by concentration of the digested peptides on Evotips to minimize losses otherwise associated with additional desalting steps. The standardized workflows can robustly process up to 192 samples in less than 10 hours with high precision and sensitivity. Close to 1,500 protein groups were quantified from neat plasma, increasing to more than 5,000 protein groups from the MagNet workflow with a throughput of 100 samples analyzed per day using the Orbitrap Astral mass spectrometer.



Figure 1: Plasma proteomics has many applications and shows great promise.

2. Method details

Plasma was extracted as per the Early Detection Research Network (EDRN) SOP³. For the neat workflow, 1 μ l plasma was diluted 180X in one-pot lysis, reduction, and alkylation buffer (1% SDS, 5 mM TCEP, 10 mM CAA in 50 mM TEAB) and subsequently 6 μ l was added to 5 μ l MagReSyn Hydroxyl magnetic beads (ReSyn Biosciences). Acetonitrile was transferred to each well of the sample plate to a final concentration of 80% driving on-bead protein capture. Post 10 min. bind step, and a single wash using 100% acetonitrile, digestion was carried out for 4 hours using a combination of 10 ng LysC and 40 ng Trypsin. Following digestion, 40% of the digest was directly loaded on Evotips on the OT-2. In the case of the Mag-Net workflow, 4 μ l plasma was mixed with 4 μ l bind buffer (100 mM Bis-Tris Propane, 150 mM NaCl pH 6.5) and 1 μ l MagReSyn SAX beads (ReSyn Biosciences), and diluted with 32 μ l wash buffer (50 mM Bis-Tris Propane, 150 mM NaCl pH 6.5). Three sequential 12 min. bind steps were followed by three washes using 100 μ l wash buffer. Next, one-pot buffer (1% SDS, 5 mM TCEP, 10 mM CAA in 50mM Tris-HCl pH 8.5) was used to solubilize, reduce, and alkylate proteins during a 1-hour on-deck incubation.

The remainder of the steps followed the neat workflow. Proteins were re-captured using acetonitrile-induced on-bead aggregation followed by a single wash and digestion for 4 hours using a combination of 75 ng LysC and 300 ng trypsin. 40% of the resulting digest was loaded on Evotip. Samples were analyzed with the 100 SPD method (EV1109 column, Evosep, operated at 40 °C) coupled to an Orbitrap Astral mass spectrometer (Thermo Scientific). Spray voltage was set to 1900 V, and the heated capillary temperature was 275 °C. The mass spectrometer was operated at a full MS resolution of 240,000 with a full scan range of 380 – 980 m/z. The full MS AGC was set to 500%. MS/MS scans were recorded with a 3 Th isolation window, 7 ms maximum ion injection time. MS/MS scanning range was 380-980 m/z. The isolated ions were fragmented using HCD with 25% NCE. Data was analyzed using DIA-NN (version 1.8.1) in library-free mode against the reviewed human proteome (Uniprot, Oct 2020, 20,600 entries) with trypsin/P as digestion enzyme allowing 2 missed cleavages. All conditions were searched separately with match between runs enabled across replicates within the same condition.

3. Maximizing efficiency

The end-to-end Mag-Net workflow was adapted from the original publication² to facilitate integrated Evotip loading and downscaling of the plasma input to reduce costs of expensive reagents such as enzymes by only digesting what is needed for a single LC-MS injection. This was implemented on the OT-2, where digestion was carried out for 4 hours at ambient temperature to ensure high throughput despite the limited deck space on the robot. 1 μ l of plasma was used as input for the neat plasma

workflow, from which ~1 μ g was digested using a PAC assisted digestion, whilst the Mag-Net workflow used 4 μ l plasma as input (Figure 2). The protocol utilizes direct loading of Evotips after digestion, which leads to increased sensitivity compared to conventional elution, dry-down, and resuspension steps. The protocol is sustainable by design as just a single pipette tip is used for all sample transfer steps and minimal reagents are used for digesting only what is analyzed.

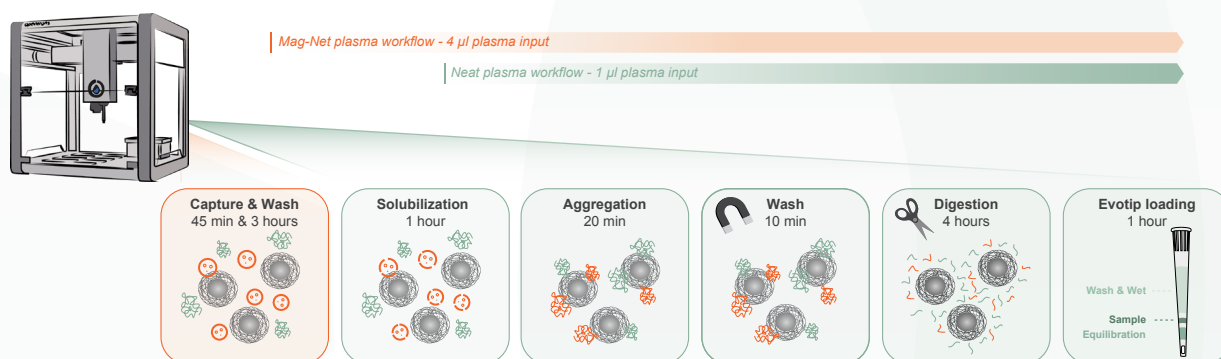


Figure 2: Schematic representation of the OT-2 workflows for neat and Mag-Net plasma profiling.

4. Reproducible workflows

We assessed the robustness of both workflows across two plates on the same OT-2, with a total of 24 samples of both neat plasma and MagNet enriched plasma, respectively (Figure 3). Using the 100 SPD method, the neat workflow consistently identified over 1,400 unique protein groups and >15,000 precursors. Importantly, 1,033 protein groups and 6,519 precursors were quantified in all samples. Using the automated Mag-Net workflow, more than 5,000 protein groups and 45,000 precursors

were quantified in each sample and 25,264 precursors and 4,060 protein groups were quantified in all samples. The precision in the data was excellent with a median coefficient of variation (CV) on protein level of approximately 10% for both neat and Mag-Net workflows. For context, the technical variability in the LC-MS setup contributed with a median CV of less than 4%, providing a solid platform for robust high-throughput analysis (Figure 3).

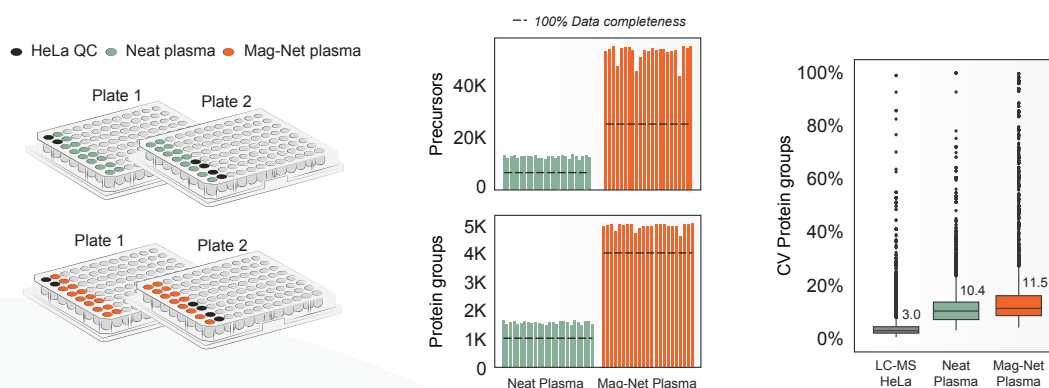


Figure 3: Experimental setup and precursor and protein ID levels. CVs representing LC-MS and workflow reproducibility. ~40% of each peptide digest was loaded on Evotips and analyzed with 100 SPD.

An asset of this strategy is the option to process both the neat and Mag-Net-based plasma in parallel from just 5 μ l of plasma. Besides the obvious proteome depth in the datasets, the two workflows are complementary from a biological perspective. The neat workflow predominantly identifies soluble, highly abundant plasma proteins. While the dynamic range of these are more than 10-fold, the sensitivity of

the method enables the identification of proteins in the ng/ml range based on estimated protein concentration levels reported in Human Disease Blood Atlas (HBDA)⁴ (Figure 4). This enables the monitoring of numerous relevant proteins including 52% (59 of 113) of FDA-approved plasma biomarkers where 72% of these were quantified with a CV<20%.

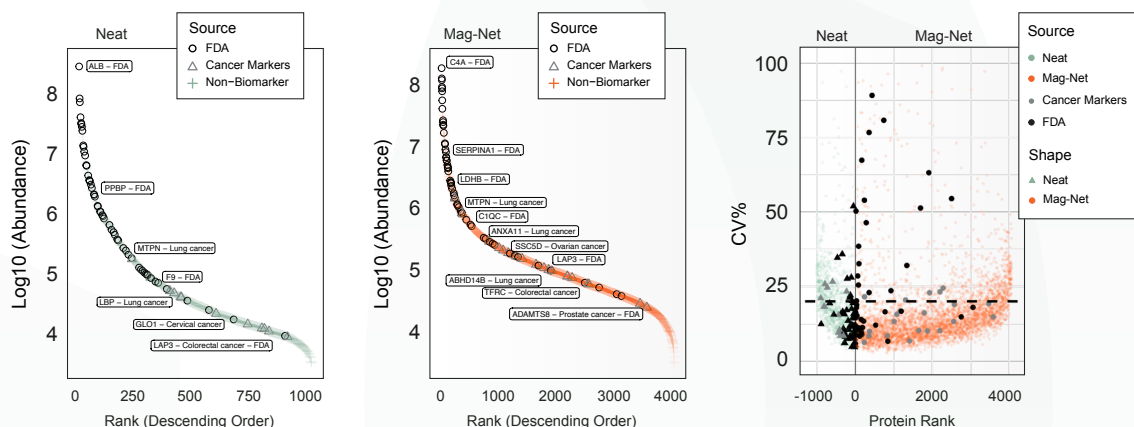


Figure 4: Neat and Mag-Net plasma dynamic range plot (descending) of 100% coverage proteins, annotated with CV% and isoform specific FDA and HBDA putative biomarkers.

The Mag-Net workflow reduces the dynamic range of the enriched sub-proteome by depleting soluble highly abundant plasma proteins whilst enriching for membrane-bound vesicles. 96 of the top 100 extracellular vesicle marker proteins listed in Vesiclepedia were identified in the Mag-Net dataset, where 82% of these were quantified with a

CV<20%. Further, numerous proteins listed as putative markers in the Human Disease Blood Atlas with estimated blood concentrations in the low pg/ml range were measured in this workflow with CVs<20%. In combination, these proteins could be utilized as reporters of nine different cancer types in relevant cancer cohorts and studies (Figure 4).

5. Conclusion

Here, we describe the development of a complete, fully automated plasma profiling pipeline that is cost-efficient by design and seamlessly integrated with high-throughput LC-MS analysis using the Evosep One. This enables the processing of 192 samples in short time - from raw plasma to ready-to-analyze Evotips. Both workflows, for neat plasma and membrane-bound vesicle enrichment provide good quantitative precision with median protein CVs of approximately 10%. These can be processed independently of each other or within one OT-2 with just 5 µl of total plasma required

for both methods. Coupling this sensitive, end-to-end analytical pipeline to the rapid 100 SPD method delivers deep plasma proteome profiles with more than 5,300 protein groups identified. This depth allowed the monitoring of low-abundant biomarkers.

The ability to tackle large clinical cohorts in a robust, cost-effective, high-throughput, and sensitive manner opens the door to next-generation protein biomarker discovery as well as expediting clinical implementation of existing marker panels by providing an efficient orthogonal assay to be used for biomarker validation.

Evosep One is for General Laboratory Use.

References

1. Batth TS., Tollenaere M., Rüther P., Gonzalez-Franquesa A., Prabhakar BS., Bekker-Jensen S., Deshmukh AS., Olsen JV (2019) Protein Aggregation Capture on Microparticles Enables Multipurpose Proteomics Sample Preparation. Mol Cell Proteomics., mcp.TIR118.001270
2. Wu CC., Tsantilas KA., Park J., Plubell D., Naicker P., Govender I., Buthelezi S., Stoychev S., Jordaan J., Merrihew G., Huang E., Parker ED., Riffle M., Hoofnagle AN., MacCoss MJ (2023) Mag-Net: Rapid enrichment of membrane-bound particles enables high coverage quantitative analysis of the plasma proteome. BioRxiv., <https://www.biorxiv.org/content/10.1101/2023.06.10.544439v1.full.pdf>
3. Tuck MK., Chan DW., Chia D., Godwin AK., Grizzle WE., Krueger KE., Rom W., Sanda M., Sorbara L., Stass S., Wang W., Brenner DE (2010) Standard Operating Procedures for Serum and Plasma Collection: Early Detection Research Network Consensus Statement Standard Operating Procedure Integration Working Group. J Proteome Res., 10.1021/pr800545q
4. Alvez MB., Edfors F., Feilitzén KV., Zwahlen M., Mardinoglu A., Edqvist PH., Sjöblom T., Lundin E., Rameika N., Enblad G., Lindman H., Höglund M., Hesselager G., Ståhlberg K., Enblad M., Simonson OE., Häggman M., Axelsson T., Åberg M., Nordlund J., Zhong W., Karlsson M., Gyllenstein., Ponten F., Fagerberg L., Uhlén M (2023) Next generation pan-cancer blood proteome profiling using proximity extension assay. Nat Commun., 10.1038/s41467-023-39765-y